

## Human Serum Pooled Donors, 3X Charcoal Stripped

<b>Catalog No.</b>	CSI20430A	<b>Quantity:</b>	100 ml
	CSI20430B		1 liter

**Description:** Charcoal is used to strip hormones and other analytes from plasma derived, defibrinated, USA origin, pooled human plasma.  
Lot-specific results reported on Certificate of Analysis for Glucose, Sodium, Potassium, Chloride, Calcium, Phosphorus, Total Protein, Albumin, Bioburden, Total Cholesterol, Triglycerides, Testosterone, pH, Cortisol, Estradiol, Total T3, Total T4, B.U.N., Creatinine, and Uric Acid.

**Appearance:** Clear yellow liquid.

**Preservative:** None

**Handling:** Thaw in a 37°C water bath. Gently mix every 10 minutes to ensure even thawing. Thawed serum should be used immediately or stored at 2-8°C for up to 24 hours. Prepare single use aliquots and store below -20°C.

**Country of Origin:** USA

**Storage & Stability:** Store at -20°C to -80°C for 1 year. Upon initial thawing, aliquot and store at -20°C to -80°C. **Avoid repeated freeze/thaw cycles.**

**Certification:** Non-reactive for HIV-1, HCV, HBV by NAT, HBsAg, HCV Ab, HIV 1 and 2 Ab, and RPR by currently approved FDA methods. However, because no test method can offer complete assurance that infectious agents are absent, this material should be handled at the Bio-safety level 2 (BSL 2) as recommended for any potentially infectious human serum or blood specimen in the CDC/NIH manual "Bio-safety in microbiological and Biometical Laboratories", 1999.

**NOT FOR HUMAN USE. FOR RESEARCH ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.**



**Cell Sciences**<sup>®</sup>  
65 Parker Street  
Unit 11  
Newburyport, MA 01950

Toll Free: 888-769-1246  
Phone: 978-572-1070  
Fax: 978-992-0298

E-mail: [info@cellsciences.com](mailto:info@cellsciences.com)  
Website: [www.cellsciences.com](http://www.cellsciences.com)